

FAQ

Ni Sepharose High Performance
HisTrap HP
HisTrap HP Kit



Q When should I use Ni Sepharose™ 6 Fast Flow instead of Ni Sepharose™ High Performance?

A Both Ni Sepharose 6 Fast Flow and Ni Sepharose High Performance are optimized for purification of histidine-tagged proteins and the media are pre-charged with Ni²⁺.

Ni Sepharose 6 Fast Flow is designed for fast, reliable scale-up of histidine-tagged protein purification. The medium is also well suited for batch/gravity flow purification and multi-well plate screening.

Ni Sepharose High Performance is optimized for high performance purification of histidine-tagged proteins with a chromatography system, such as ÄKTAdesign™ from GE Healthcare.

Q What is the protein binding capacity of Ni Sepharose High Performance?

A *Dynamic binding capacity:* At least 40 mg (histidine)₆-tagged protein/ml medium

Dynamic binding capacity conditions:

Sample: 1 mg/ml (histidine)₆-tagged pure proteins (M_r 28 000 or M_r 43 000) in binding buffer (QB_{10%} determination) or (histidine)₆-tagged protein bound from *E. coli* extract

Column volume: 0.25 ml or 1 ml

Flow rate: 0.25 ml/min or 1 ml/min

Binding buffer: 20 mM sodium phosphate, 0.5 M NaCl, 5 mM imidazole, pH 7.4

Elution buffer: 20 mM sodium phosphate, 0.5 M NaCl, 500 mM imidazole, pH 7.4

Note: Dynamic binding capacity is protein dependent.

Q What are the general running conditions recommended?

A Recommended starting conditions:

Binding buffer: 20 mM sodium phosphate, 0.5 M NaCl, 20–40 mM imidazole, pH 7.4

Elution buffer: 20 mM sodium phosphate, 0.5 M NaCl, 500 mM imidazole, pH 7.4

Note: The concentration of imidazole that will give optimal purification results (in terms of purity and yield) is protein-dependent, and is usually slightly higher for Ni Sepharose High Performance than for similar IMAC media on the market. This can be useful during the washing step as a higher concentration of imidazole in the wash buffer may wash out contaminants that have bound to the medium. This can result in elution of a more pure target protein.

Higher concentration of imidazole will not cause any problems in terms of increased baseline absorbance if a high purity imidazole is used, such imidazole gives essentially no absorbance at 280 nm.

Q Will the high capacity of Ni Sepharose High Performance result in higher non-specific binding?

A The non-specific binding of proteins is a result of several factors. Other proteins will only bind more to Ni Sepharose High Performance if these non-specific proteins are interacting directly with the Ni²⁺ or the metal chelating ligand. In addition, non-specific proteins can potentially be washed off at higher concentrations of imidazole, while keeping the target protein bound to the medium.

Q With the higher binding capacity, do you have to use harsh conditions to remove the target protein?

A Harsh conditions are not necessary to remove most target proteins. The proteins tested are all eluted below 500 mM imidazole; frequently between 100-300 mM imidazole. The protein is efficiently eluted from the column, resulting in high yields.

Q How many times can you run the Ni Sepharose High Performance while maintaining the same capacity?

A This is sample dependent. Reusing Ni Sepharose High Performance depends on the nature of the sample. To prevent cross-contamination, only reuse the column when purifying identical histidine-tagged proteins.

Q How do you measure the amount of Ni²⁺ in the eluent?

A This quantitative measurement assay is as follows:

The medium is run with 10 column volumes of a buffer with pH 4.0 and then the Ni²⁺ still bound on the column is measured spectroscopically and compared with the amount Ni²⁺ loaded.

Q Can you strip and regenerate the Ni Sepharose High Performance?

A Ni Sepharose High Performance can be stripped with 50 mM EDTA in 20 mM sodium phosphate, 0.5 M NaCl, pH 7.4 and recharged with Ni²⁺ or other divalent cations, such as Cu²⁺, Co²⁺, Zn²⁺, etc.

Q How can I strip and recharge Ni Sepharose High Performance?

A Recommended stripping buffer: 20 mM sodium phosphate, 0.5 M NaCl, 50 mM EDTA, pH 7.4.

1. Strip the column by washing with 5 to 10 column volumes of stripping buffer.
2. Washing with 5 to 10 column volumes of binding buffer.
3. Wash with 5 to 10 column volumes of distilled water.
4. Re-charge the column.

To recharge the water-washed column:

1. Load 0.5 column volumes of 0.1 M metal salt solution in distilled water on the column. Metal chloride and sulfate salts (e.g., 0.1 M NiSO₄) are commonly used.
2. Wash with 5 column volumes of distilled water.
3. Wash with 5 column volumes of binding buffer.

Note: It is important to wash with binding buffer as the last step to get the correct pH.

Q How is the Ni²⁺ leakage at low pH?

A At pH 4, the loss of Ni²⁺ is < 5% on Ni Sepharose High Performance (reproducible) compared to average 9% for Ni-NTA Superflow from Qiagen (wide batch to batch variations).

Note: Loss of metal ions is more pronounced at lower pH.

Q My target protein is extremely sensitive to Ni²⁺. How do I reduce the Ni²⁺ leakage to a real minimum?

A Leakage of Ni²⁺ from Ni Sepharose High Performance is low under all conditions and resistance towards reducing agents is thus high.

Note: If using buffers/sample including reducing agents it is recommended to perform a blank run without reducing agents in the buffers before loading sample (to remove any weakly bound Ni²⁺ ions):

Blank run: Use binding and elution buffer *without* reducing agents

1. Wash the column with 5 column volumes of distilled water.
1. Wash with 5 column volumes of binding buffer.
2. Wash with 5 column volumes of elution buffer.
3. Equilibrate with 10 column volumes of binding buffer.

Q What different additives can I use together with Ni Sepharose High Performance and HisTrap™ HP?

A Tests performed in our laboratories show that Ni Sepharose High Performance is compatible with the following compounds at the concentrations given.

Reducing agents	5 mM DTE 5 mM DTT 20 mM β-mercaptoethanol 5 mM TCEP 10 mM reduced glutathione	
Denaturing agents	8 M urea* 6 M guanidine hydrochloride*	
Detergents	2% Triton™ X-100 2% Tween™ 20 2% NP-40 2% cholate 1% CHAPS	Non-ionic detergent Non-ionic detergent Non-ionic detergent Anionic Detergent Zwitterionic Detergent
Additives	500 mM imidazole 20% ethanol 50% glycerol 100 mM Na ₂ SO ₄ 1.5 M NaCl 1 mM EDTA** 60 mM citrate**	
Buffer solutions	50 mM sodium phosphate, pH 7.4 100 mM Tris-HCl, pH 7.4 100 mM Tris-acetate, pH 7.4 100 mM HEPES, pH 7.4 100 mM MOPS, pH 7.4 100 mM sodium acetate, pH 4*	

* Tested for 1 week at +40 °C

** Generally, chelating agent should be used with caution (and only in the sample, not in buffers). Any metal ion stripping may be counteracted by addition of a small excess of MgCl₂ before centrifugation or filtration of the sample. Note that stripping effects may vary with applied sample volume.

Q How do I remove imidazole from my target protein after elution?

A If imidazole needs to be removed from the protein, use HiTrap™ Desalting 5 ml (can be coupled in series to increase the sample volume), a PD-10 column, or HiPrep™ 26/10 Desalting, depending on the sample volume.

Q Do you see an improved purity (less non-specific binding) with high concentrations of NaCl?

A In most cases, concentration of ions, imidazole, and other additives in the binding and wash buffers will influence the level of non-specific binding. Higher concentrations of NaCl (500 mM) generally reduce the level of non-specific binding.

Q Is it possible to use high concentrations of Tris-HCl buffer and how does this compatibility to Tris buffers compare to other suppliers?

A These are the Tris buffers systems that have been tested and work well.

50 mM Tris-HCl, pH 7.4

100 mM Tris-HCl, pH 7.4

100 mM Tris-acetate, pH 7.4

Tris buffers in high concentrations can adversely affect the binding of histidine-tagged target proteins to nickel medium from other suppliers; therefore, lower concentrations of Tris buffers (less than 50 mM) or phosphate buffers are generally used.

Q How can I clean Ni Sepharose High Performance?

A Note: Strip off the metal ions before cleaning (to prevent precipitation of metal salt)

To remove precipitated proteins, hydrophobically bound proteins and lipoproteins:

1. Washing the column with 1 M NaOH, contact time usually 1 to 2 hours (12 hours or more for endotoxin removal).
2. Wash the column with approximately 10 column volumes of binding buffer.
3. Wash with 5 to 10 column volumes of distilled water.

To remove strongly hydrophobically bound proteins, lipoproteins and lipids:

1. Wash the column with 5 to 10 column volumes 30% isopropanol for about 15 to 20 minutes.
2. Wash the column with approximately 10 column volumes of distilled water.
3. Re-charge the medium or wash with 5 column volumes 20% ethanol for storage.

Alternatively:

1. Wash the column with 2 column volumes of detergent in a basic or acidic solution. Use, for example, 0.1% to 0.5% nonionic detergent in 0.1 M acetic acid, contact time 1 to 2 hours.
2. After treatment, always remove residual detergent by washing with 5 column volumes of 70% ethanol.
3. Wash the column with approximately 10 column volumes of distilled water.
4. Re-charge the medium or wash with 5 column volumes 20% ethanol for storage.

To remove ionically bound proteins:

1. Wash the column with 1.5 M NaCl solution, contact time 10 to 15 minutes.
2. Wash the column with approximately 10 column volumes of distilled water.
3. Re-charge the medium or wash with 5 column volumes 20% ethanol for storage.

Q Is it possible to use Ni Sepharose High Performance by gravity or in 96- or 384-well multiwell plates for screening?

A Yes it is possible, but for gravity usage we recommend using Ni Sepharose 6 Fast Flow.

www.gehealthcare.com/his

GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala
Sweden

ÄKTAdesign, HiPrep, HisTrap, and Sepharose are trademarks of GE Healthcare companies. GE, imagination at work, and GE monogram are trademarks of General Electric Company.

Triton is a trademark of Union Carbide Chemicals and Plastics Co. Tween is a trademark of ICI Americas Inc. Qiagen is a trademark of the Qiagen group. Superflow is a trademark of Sterogene Bioseparations Inc.

All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare that supplies them. General Electric Company reserves the right, subject to any regulatory and contractual approval, if required, to make changes in specifications and features shown herein, or discontinue the product described at any time without notice or obligation. Contact your local GE Healthcare representative for the most current information.

© 2006 General Electric Company – All rights reserved.

GE Healthcare Bio-Sciences AB, a General Electric Company.

GE Healthcare Bio-Sciences AB
Björkgatan 30, 751 84 Uppsala, Sweden

GE Healthcare Europe GmbH
Munzinger Strasse 5, D-79111 Freiburg, Germany

GE Healthcare UK Ltd
Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK

GE Healthcare Bio-Sciences Corp
800 Centennial Avenue, P.O. Box 1327, Piscataway,
NJ 08855-1327, USA

GE Healthcare Bio-Sciences KK
Sanken Bldg. 3-25-1, Hyakunincho, Shinjuku-ku,
Tokyo 169-0073, Japan

Asia Pacific T +85 65 62751830 F +85 65 62751829 • Australasia T +61 2 8820 8299 F +61 2 8820 8200 • Austria T 01 /57606 1613 F 01 /57606 1614 • Belgium T 0800 73 890 F 02 416 8206 • Canada T 1 800 463 5800 F 1 800 567 1008 • Central & East Europe T +43 1 972 720 F +43 1 972 722 750 • Denmark T +45 70 25 24 50 F +45 45 16 2424 • Eire T 1 800 709992 F +44 1494 542010 • Finland & Baltics T +358 9 512 3940 F +358 9 512 39439 • France T 01 69 35 67 00 F 01 69 41 98 77 • Germany T 0800 9080 711 F 0800 9080 712 • Greater China T +852 2100 6300 F +852 2100 6338 • Italy T 02 26001 320 F 02 26001 399 • Japan T 81 3 5331 9336 F 81 3 5331 9370 • Korea T 82 2 6201 3700 F 82 2 6201 3803 • Latin America T +55 11 3933 7300 F +55 11 3933 7304 • Middle East & Africa T +30 210 96 00 687 F +30 210 96 00 693 • Netherlands T 0800-82 82 82 1 F 0800-82 82 82 4 • Norway T +47 815 65 777 F +47 815 65 666 • Portugal T 21 417 7035 F 21 417 3184 • Russia, CIS & NIS T +7 495 956 5177 F +7 495 956 5176 • Spain T 902 11 72 65 F 935 94 49 65 • Sweden T 018 612 1900 F 018 612 1910 • Switzerland T 0848 8028 10 F 0848 8028 11 • UK T 0800 515 313 F 0800 616 927 • USA T +1 800 526 3593 F +1 877 295 8102

