



## Protocol for Total RNA Purification from Fresh Cultured CHO (Chinese Hamster Ovary) Cells Using Pall Nucleic<sup>®</sup> Acid Binding Nanosep<sup>®</sup> Centrifugal Device

### 1. Consumables and Reagents

**Table 1**

*Consumables for Total RNA Purification (nuclease-free consumables are recommended)*

Supplier	Product Description	VWR Cat. No.
Pall Laboratory	Nucleic Acid Binding (NAB) Nanosep Centrifugal Device	76360-454 , 76360-456
VWR	Ethanol (not denatured)	71001-866
VWR	Spectrophotometer Cuvettes ~100 µL (260/280 nm)	47743-840
VWR	Tubes 15 mL (RNase-DNase free)	89401-574
VWR	Tubes 50 mL (RNase-DNase free)	89401-572
VWR	Microcentrifuge Tubes 1.5 mL (RNase-DNase free)	76005-210
VWR	Needles (20) (0.9 mm)	89219-312
VWR	Syringe (10 mL)	89215-218

**Table 2**

*Reagents for Total RNA Purification (nuclease-free reagents are recommended)*

Supplier	Product Description	VWR Cat. No.
Various	Proteinase K	N/A
Various	DNaseI set	N/A
Various	Lysis Buffer	N/A
Various	Wash Buffer 1	N/A
Various	Wash Buffer 2	N/A
Various	Nuclease-free water	N/A
Various	TE Buffer (pH 7.5)	N/A
VWR	Tris Buffer pH 7.0 (1 M)	89500-584
VWR	Lysozyme	97062-136
VWR	DTT (25 g)	97063-758

## 2. Instruments

- Microcentrifuge
- Spectrophotometer
- Vortex

## 3. Important Points Before Starting

- Clean all equipment/material to be used for RNA extraction.
- All centrifugation steps are performed at room temperature at 10,000 – 14,000 x g.
- It is essential to work quickly and efficiently when working with RNA.
- For each NAB Nanosep device insert there are three receiver tubes. This is enough to complete the below process. Use only the supplied receiver tubes with the NAB Nanosep device.
- All buffers should be allowed to equilibrate to room temperature before use.
- Briefly centrifuge tubes after vortexing to remove drops from inside the lid.
- Change pipette tips between all liquid transfers. Pall recommends use of sterile RNA-free pipette tips.

## 4. Protocol

1. Estimate the quantity of cells before starting. CHO cell concentrations up to  $10^7$  may be used.
2. Centrifuge the culture at 1,000 x g for 10 minutes to pellet the appropriate number of cells. Remove the supernatant carefully by aspiration.
3. Add 500  $\mu$ L of lysis buffer (supplemented with DTT) per tube. Vortex the mixture vigorously.
4. Using a blunt 20-gauge needle fitted to an RNase-free syringe, pass the lysate in and out of the syringe ~10 times followed by 20 seconds of vigorous vortexing.
5. Add 500  $\mu$ L of 100% non-denatured ethanol to the homogenised lysate. Mix well by pipetting but DO NOT centrifuge.
6. Transfer up to 500  $\mu$ L of the lysed cells to the NAB Nanosep device insert inside a receiver tube, including any precipitate. Close the lid and centrifuge for 60 seconds at 10,000 – 14,000 x g. Discard the flow-through but re-use the collection tube for the next step.
7. Repeat the previous step if samples are greater than 500  $\mu$ L using the same device.
8. Optional DNase digestion steps: (If you do not require DNase digestion step move directly to step 9) (See supplier instructions for DNase I preparation)
  - a. Add 350  $\mu$ L of wash buffer 1 to the NAB Nanosep device insert. Close the lid and centrifuge for 60 seconds at 10,000 – 14,000 x g to wash the membrane.
  - b. Discard the flow-through and retain the receiver tube for the next step.
  - c. Add 80  $\mu$ L of the prepared DNase I solution directly on to the NAB Nanosep device filter membrane.
  - d. Incubate at room temperature for at least 15 minutes.
  - e. Add 350  $\mu$ L wash buffer 1 to the NAB Nanosep device. Close the lid and centrifuge at 10,000 – 14,000 x g for 60 seconds. Discard the flow-through, retain the receiver tube and proceed to step 10.

9. Wash the NAB Nanosep device membrane with 500  $\mu$ L of wash buffer 1 and centrifuge for 1 minute at 10,000 – 14,000 x g. Discard the flow through and retain the receiver tube for the next step.
10. Add 500  $\mu$ L of wash buffer 2 to the NAB Nanosep device, close the lid and centrifuge for 60 seconds at 10,000 – 14,000 x g. Ensure all the solution has passed through the filter membrane to avoid any carryover.
11. Repeat step 10 but centrifuge for 2 minutes.
12. Carefully remove the NAB Nanosep device insert, being careful not to allow the filtrate to contact the insert, and discard the receiver tube.
13. Place the NAB Nanosep device insert into a clean receiver tube (provided), close the lid and centrifuge for 60 seconds at 10,000 – 14,000 x g.
14. Discard the filtrate tube and place the NAB Nanosep device insert into a clean receiver tube.
15. Add 50  $\mu$ L of RNase-free water directly on to the NAB Nanosep device insert filter membrane. Close the lid and incubate at room temperature for 1 minute. Centrifuge for 60 seconds at 10,000 – 14,000 x g to elute the RNA from the filter membrane.
16. Optional: Repeat the elution step with a further 50  $\mu$ L RNase-free water in the same device, in the same receiver tube.

### Storage of RNA

Purified RNA can be stored in RNase-free water at  $-20\text{ }^{\circ}\text{C}$  or  $-70\text{ }^{\circ}\text{C}$  for 1 year.

### Quantification of RNA

RNA concentration can be determined by measuring the absorbance at 260 nm ( $A_{260}$ ) in a spectrophotometer (see details below). For small quantities of RNA however, it can be difficult to determine these amounts photometrically. Smaller quantities of DNA can be accurately quantified using fluorometric quantification.

### Spectrophotometric quantification of RNA

$A_{260}$  readings should be greater than 0.15 to ensure significance. An absorbance reading of 1.0 at 260 nm corresponds to 44  $\mu$ g of RNA per mL. This is only valid for measurements at neutral pH however. As a result, if it is necessary to dilute the RNA sample, ensure that the dilution buffer is of neutral pH.



Ordering:  
vwr.com

© 2020 Pall Corporation. Pall (PALL) and Nanosep are trademarks of Pall Corporation. ® indicates a trademark registered in the USA.  
*Filtration. Separation. Solution.* is a service mark of Pall Corporation.

4/20, PDF, GN19.1112

*Filtration. Separation. Solution.*<sup>SM</sup>

800 932 5000 | VWR.COM

Prices, product, and/or services details are current when published and subject to change without notice. | Certain products or services may be limited by federal, state, provincial, or local regulations. | VWR, part of Avantor, makes no claims or warranties concerning sustainable/green products. Any claims concerning sustainable/green products are the sole claims of the manufacturer and not those of VWR International, LLC and/or Avantor, Inc. or affiliates. All prices are in US dollars unless otherwise noted. Offers valid in US and Canada, void where prohibited by law or company policy, while supplies last. | Trademarks are owned by Avantor, Inc. or its affiliates, unless otherwise noted. | Visit vwr.com to view our privacy policy, trademark owners, and additional disclaimers. © 2020 Avantor, Inc. All rights reserved.

0320 Lit. No. 160144W